


## Home Screen Reset

Technical Note 163

## Introduction

The DeNovix® DS-11 software is an Android OS based system with a custom home screen. On rare occasions, the Home screen may not be properly displayed after a software update.

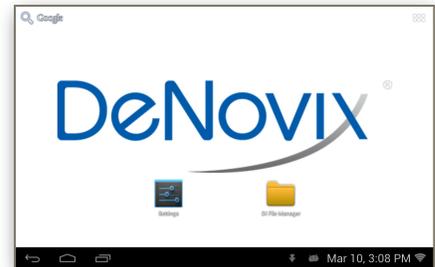


Incorrect Home Screen

## Procedure

Follow the steps below to reset the software to the proper Home screen:

1. Tap the 6 dot **More** icon found at the top right of the screen
2. The Android formatted application screen will appear. Swipe to the second page and launch the **Settings** app.
3. Scroll down and select **Apps**.
4. Scroll down and select **Launcher** from the list.
5. Scroll down and select **Clear defaults**.
6. Return to the Home screen using the bottom navigation bar **Home** icon indicated by the circle in the image to the right.
7. Complete action by selecting **DeNovix DS11 Home** and **Always**.

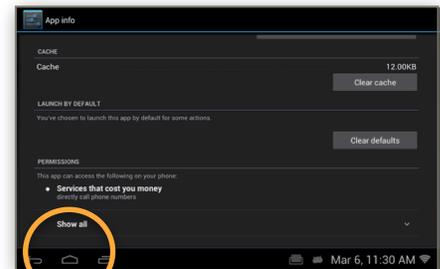

Home Screen Icon

Please contact Customer Support at [info@denovix.com](mailto:info@denovix.com) or 302-442-6911 for additional information. Outside of the US, please contact your local distributor.