


## Introduction

The software for the DeNovix® Spectrophotometer / Fluorometer Series instruments enables printing of data and screen captures via network printers. This document provides a step-by-step guide to exporting from a DeNovix instrument to a network printer.

## Procedure for Network Printer Setup

1. Download and install the most current version of DeNovix software to ensure that the Network Printing app includes the most recent drivers.
  - Note for Ethernet users: Ensure that Wi-Fi is turned off in the Settings app.
2. Using any measurement app or the Data app, select the data of interest and use the top right drop down menu to export data or a screen capture.
3. Select **Network Printer** from the export option listed.
4. Tap **Print** or **Printer Not Selected**, depending on which of the configurations below is shown. If a pop up asking to “Setup printer now” appears, tap **OK**.

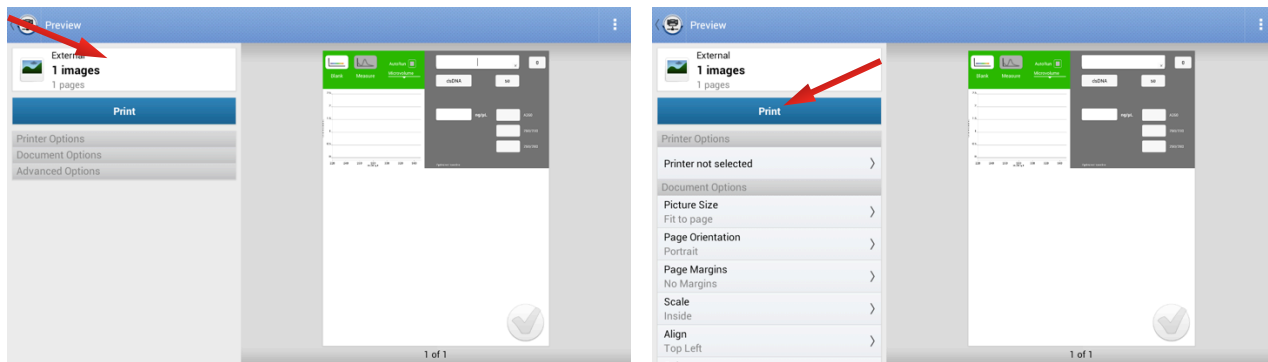


5. The Manage Printers screen should appear as below. Tap **Network Printers**.




## Procedure for Network Printer Setup continued

- The available printers will be displayed as shown below. Search for and select the printer of interest.
  - Note: The Scan for Wi-Fi Printers button searches for all printers on the network, including those that are ethernet-connected.

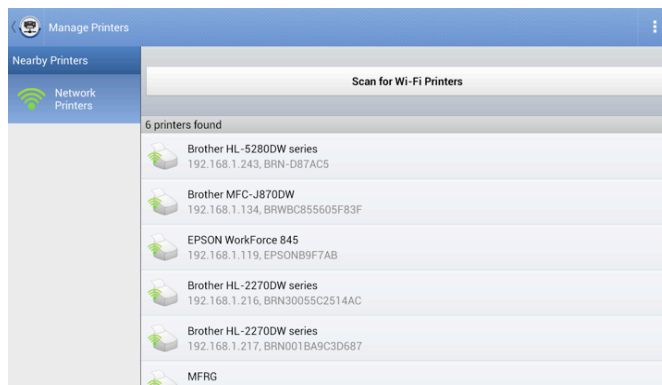

- Select the appropriate print settings using the options list on the left of the screen.
- Tap **Print** to print the selection.

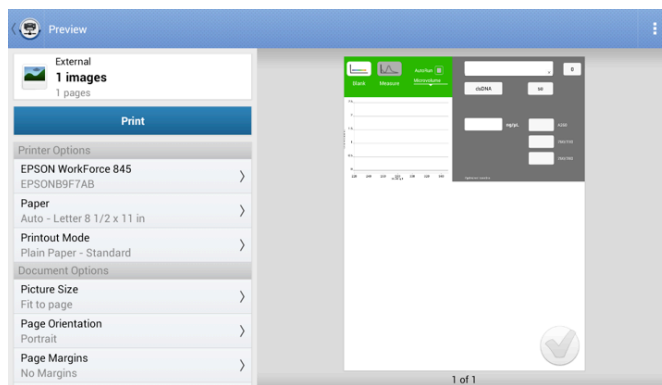

The Network Printing feature is compatible with 90% of commercial printers. Below is a list of printers that are known to work successfully with DeNovix instruments.

- Brother HL-2270DW
- Brother HL-4570CDW
- Brother HL-5280DW
- Epson L355
- Epson 845
- HP Color LaserJet CM1312nfi MFP
- HP Color LaserJet 4500
- HP Deskjet 3050A
- HP Officejet 6600
- HP Officejet 6700 Premium

Please contact Customer Support at [info@denovix.com](mailto:info@denovix.com) or 302-442-6911 for additional information. Outside of the US, please contact your local distributor.

Please see the DS-11 /FX Series Troubleshooting Guide available at [www.denovix.com](http://www.denovix.com) for troubleshooting guidance.

A DYMO® USB label printer is a space-saving, direct-connected alternative to a networked printer.

*DYMO® is a registered trademark of DYMO Corporation.*